

GenePasQ Data Sheet / Reading Guide
(AHPND/EMS, IHHNV, WSSV, EHP)

Title: _____

Date: ()

Name: ()

• Gene Amplification by PCR

Multiplex PCR Plus Kit


(Lot No.) _____

①	95 °C	5 min
②	95 °C	15 sec
③	60 °C	30 sec
④	72 °C	15 sec
⑤	4 °C	


35 cycles

• GenePasQ

(Lot No.) _____

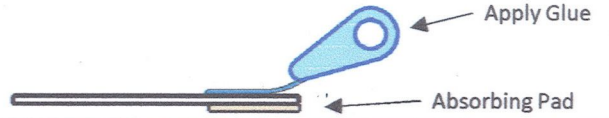

Paste the GenePasQ Strips by applying glue behind the Absorbing Pad of the Strips.

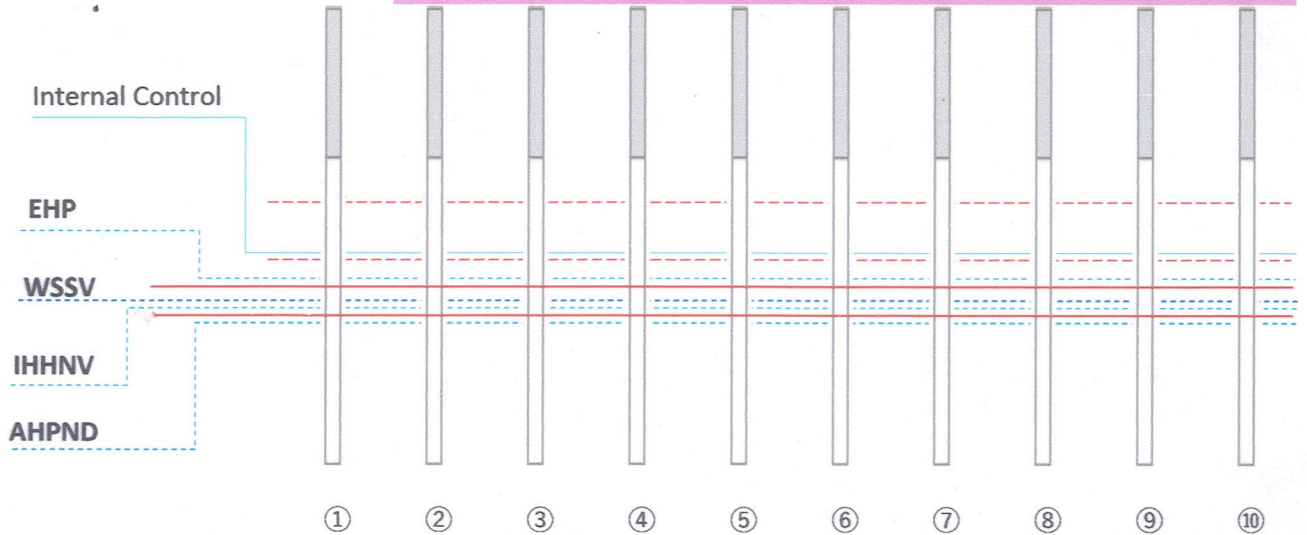

EHP										
WSSV										
IHHNV										
AHPND										

** Internal control (IC) line often becomes faint when a target gene is amplified.*

Test		Comments
①		
②		
③		
④		
⑤		
⑥		
⑦		
⑧		
⑨		